

ISO - Flex

***The Flexible
Stimulus Isolator***



optically isolated

***constant current /
constant voltage***

bipolar pulses

5-Year Warranty

ISO - Flex


The Flexible Stimulus Isolator



Flexibility

ISO-Flex is a flexible stimulus isolator since:

- ISO-Flex is optically-isolated.
- You can switch the unit to constant current or constant voltage using the CURRENT/VOLTAGE switch.
- You get bipolar pulses by connecting two isolators. The connection is as follows:

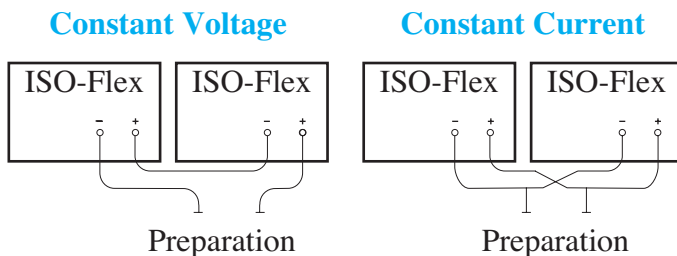

Each isolator should receive its input at its proper time. One input to drive the positive phase of the bipolar pulse, and the other to drive the negative phase.

In the constant voltage configuration both isolators should be set to constant voltage and in the constant current configuration both isolators should be set to constant current.

ISO-Flex is supplied with 2 batteries: a standard 9V battery in the ISO-Flex box and a 90V battery in the battery housing.

The 9V battery is used for the operation of the isolator. The 90V battery is used to drive the output pulses.

Set Up

To set up the ISO-Flex connect its voltage wire to the battery housing.

To reduce stimulus artifact and other noise:

Use short wires for the stimulus outputs (locate the ISO-Flex unit close to your preparation). Twist the wires together.

It is recommended that the wires not be shielded in order to eliminate capacitance to ground.

The metal chassis of the battery housing shields the 90V battery. Ground it via its banana jack.

Specifications

Input: 5-10V

Voltage output: 0-90V

Current output: 0-10mA

The output is divided into 3 ranges.

Regulatory compliance: UL and CE

ISO-Flex has not been approved for direct patient use.

5-Year Warranty



A.M.P.I., 123 Uziel St. P.O.B. 16477
Jerusalem 91163, ISRAEL
PHONE: +972 2 643-9338
FAX: +972 2 643-1833
E-mail: info@ampi.co.il
<http://www.ampi.co.il>